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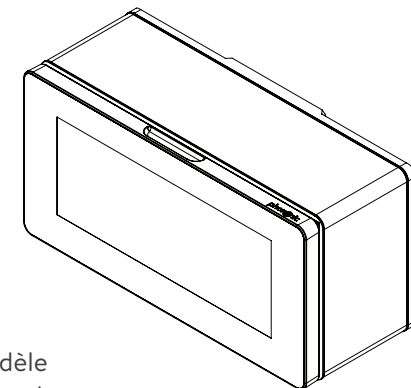
Phononic C200/F200 Series Refrigerators & Freezers

Quick Start Guide

This Quick-Start Guide applies to the following Phononic models:

C200 Merchandising Refrigerators (solid or glass door)

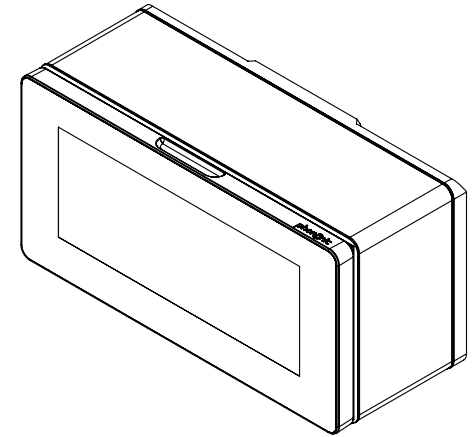
F200 Merchandising Freezers (solid or glass door)

C200-MB Laboratory & Medical Micro Benchtop Refrigerators

F200-MB Laboratory & Medical Micro Benchtop Freezers

For detailed instructions and safety precautions for your specific model, please refer to the user manual available for download at

Phononic.com/support/documentation

Step 1: Unboxing and Contents

- Move the box near install location before unpacking. *It is recommended that two people remove the unit from the packaging.*
- Inspect your unit for any shipping damage. Do not plug in the unit if there are any signs of damage.
- Contents should include the refrigerator or freezer unit, one power supply and clear storage bin (installed).
- If there are missing contents or damage to your unit, contact Phononic Customer Care: **Support@Phononic.com** or **+1.844.476.4202**

Step 2: Placement and Set Up

- Place the unit on a clean, flat, indoor surface.
- Do not block or cover the ventilation on the back of the unit.
- For best performance, we recommend a minimum rear clearance of 45cm for the F200 freezers and 10cm for the C200 refrigerators. For top/side clearance, we recommend 5cm for all models.

If your unit has not been pre-installed with decorative magnets, follow these application steps:

- Center the long (top/sides) magnet on top and wrap over the sides.
- Turn the unit over, remove the 4 feet (bumper or suction cups).
- Remove the rails from the unit with a screwdriver. Place the bolt and washer aside.
- Align magnet holes with the holes on the bottom of unit.
- Place washer on the hole and screw in the rails.
- Re-install the feet and turn unit back over.
- Apply front door magnet (if included, solid door only).

Step 3: Power On and Operation

- Connect the unit power supply cord to a 110V AC wall outlet.
- Once plugged in, the unit will automatically begin cooling.
- **There is no on-off power switch.**
- Do not load product until the default set-point temperature is reached.
- Once desired set-point temperature is reached, product may be loaded directly into the bin.
- The table shows approximate time to reach the pre-programmed set-point temperature in an ambient temperature of 24°C / 75°F.

Product	Programmed Set Temp	Approximate Pull Down Time
C200	3.3° C	2 hours
C200-MB	2.5° C or 4° C (model dependant)	2 hours
F200	-18° C	6 hours
F200-MB	-20° C	6 hours